

Thank you for your interest in our single cell analysis service. To establish a fluent cooperation please note the following points.

In case of any questions please do not hesitate to contact us.

Your Advalytix Service Team

In advance

To allow the service to take place without delay please announce your order 2-3 days in advance.

Polar body / single cell deposition

- Tear open aluminium pouch of the single cell slide..
- Take out the slide without touching the surface
- Place slide on a clean and dust free surface. Ensure that the engraved markings are readable
- Please refer to the document „ Example protocol polar body deposition“ for further information.

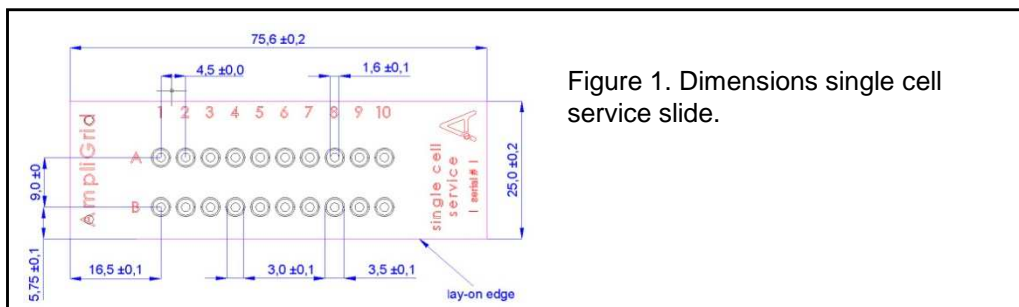


Figure 1. Dimensions single cell service slide.

- The single cell slide allows for the deposition of 20 cells max. To avoid mixing up of samples please always note the exact position of the cells and the serial number of the slide
- Cells deposited outside the engraved markings can not be analyzed!

Shipping information on the back

Shipping

- Pack the box with the single cell slide in a A4 envelope.
- Place the filled order form in the same envelope.
- Shipment of the slide with the forwarder agreed with Advalytix on the clients expense.
- Arrival at Munich central station or airport until 18:30.
- Transfer of the slide organized by Advalytix
- Please do immediately call or fax the exact arrival time to Advalytix .
Fax +49-(0)89-628366-10. Please send both pages of the order form.

Analysis

- Analysis of the cells is carried out at Advalytix.
- Given that the slide arrives until 18:30 the results will be transmitted until 9:00 the next day.
- Results are transmitted by fax followed by a hardcopy.
- Example results:

| Pos. A5 | | |
|-----------|---------|--------|
| Chromosom | present | absent |
| 1 | ✓ | |
| 2 | | ✓ |
| 3 | ✓ | |
| . | | |
| . | | |
| . | | |
| 22 | ✓ | |
| X | ✓ | |